

# 1 Heated circulating baths

## **Optima™ range**

Heated circulating baths

## **HED series**

High temperature baths/circulators

## **Gerber bath**

For milk testing and deeper tank requirements



## Heated circulating baths » Optima™ range

# Optima™ heated circulating baths and circulators

A cost-effective range of multi-purpose systems combining Grant's legendary quality and reliability. Precise temperature control for a wide range of laboratory applications.

- **Accurate and safe temperature control** – for samples and users
- **Intuitive programming and thoughtful design features** – makes working with Grant heated circulating baths and circulators easy
- **Robust, durable construction** – for longevity, reliability and long-term low cost of ownership
- **A complete range** – 32 models to cover basic through to sophisticated needs, each model represents excellent value for money



## Model selection (operating temperature)

Any of the four Grant Optima™ digital heating circulators can be combined with any of eight Grant tanks (five stainless steel and three plastic) to provide a choice of 32 models. The colour-coded summary table on page 1.6 shows you the temperature range of each combination.

The following pages showcase examples of popular combinations for different requirements.

### Liquids

We recommend the following liquids for use in Grant baths:

- 30°C to 30°C: 50% water 50% antifreeze (inhibited ethylene glycol)
- 0°C to 30°C: 80% water 20% antifreeze (inhibited ethylene glycol)
- 5°C to 99.9°C: Water
- 70°C to 150°C: Silicone fluid (viscosity ~20cS, flash point  $\geq 230^{\circ}\text{C}$ , fire point  $\geq 280^{\circ}\text{C}$ )
- 70°C to 200°C: Silicone fluid (viscosity 50cS centistokes, flash point  $\geq 285^{\circ}\text{C}$ , fire point  $\geq 340^{\circ}\text{C}$ )

## Heated circulating baths » T100, TC120, TX150 and TXF200

# Heating circulators

## T100, TC120, TX150, TXF200

The versatile Optima™ heating circulator range consists of 4 models - two general purpose: T100 and TC120 and two advanced models: TX150 and TXF200. Combine any of the four models with a Grant stainless steel or plastic tank or use independently with a clamp.

General purpose digital		Advanced digital	
T100 ambient +5 to 100°C*	TC120 ambient +5 to 120°C*	TX150 ambient +5 to 150°C*	TXF200 ambient +5 to 200°C*



T100 / TC120		TX150 / TXF200	
Features	Benefits	Features	Benefits
Stability $\pm 0.05^{\circ}\text{C}$	Excellent temperature stability and temperature control for demanding applications	Stability $\pm 0.01^{\circ}\text{C}$	Excellent temperature stability and temperature control for demanding applications
Clear, bright 4 digit LED display	Easy to view from a distance for instant reassurance of unit status	Large, bright full colour display	All key parameters visible on home screen for instant reassurance of unit status
Simple, intuitive user interface: dial and two function buttons	Easy and quick to set temperature and access menus. Minimal product training required	Icon driven home screen via a dial and two function buttons	Intuitive, quick and easy, language independent
Integral pump for external circulation (TC120)	Circulation of temperature control fluids to external apparatus / equipment	High performance integral pump for external circulation. TXF200 has variable speed	Conveniently circulate temperature control fluids to external apparatus / equipment
Model available with/without clamp (T-clamp)	Conveniently converts vessels into stirred bath, offering excellent versatility	Programming/temperature profiling (TX150, 1 program with 30 segments, TXF200 10 programs with 100 segments)	Easy and quick to configure temperature profiles to suit basic and advanced applications. Programming direct on TXF200
Low-liquid detection (float switch)	Unit will cut-out when liquid level is too low for operation	Model available with/without clamp (T-clamp)	Conveniently converts vessels into stirred bath, offering excellent versatility
User adjustable over temperature dial (TC120)	Independent safety feature and sample protection	Low-liquid detection (float switch)	Unit will cut-out when liquid level is too low for operation. Peace of mind that the unit will safely operate unattended
Fixed over temperature (T100)	Independent safety feature	5 point user calibration	Calibrate the TX150/TXF200 at any 5 temperatures against a precision reference thermometer. Provides optimum accuracy at temperatures important to the user.
Visual alarm	Alerts you when your attention is required	User adjustable over temperature dial	Independent safety feature and sample protection
2 point user calibration	Provides optimum accuracy at temperatures important to the user	Display with a choice of 5 languages (EN, DE, FR, ES & IT)	
Countdown timer (TC120)	Offers convenient reaction timing	USB/RS232 interface	Allows connection to PC or laptop for programming or data logging

- Applications:**
- Clinical, Microbiology and Pathology labs - media tempering, thawing & incubating samples
  - University research - temperature control of spectrophotometers & refractometers and jacketed vessels
  - Industrial labs - temperature probe calibration, water analysis, QC testing product, petrochemical testing, material testing, milk sample testing

Heated circulating baths » TC120-ST12 mid range showcase

showcase 1 – mid range example

Model TC120-ST12\* range 0°C to 120°C\*\*, stability ±0.05°C

Versatile mid-range model with digital thermostatic control unit and stainless steel tank and a comprehensive specification to suit most applications for precision temperature control.

- Optima™ digital thermostat (TC120) for precise temperature control
- Integral pump for external fluid circulation
- Cooling/heating range 0°C to 120°C\*\*
- Stability ±0.05°C
- 3 programmable temperature presets
- Easy to use rotary dial and two function keys

TC120-ST12 model shown

- Countdown timer with audible alarm – alerts you when your attention is required
- Simple-to-use rotary dial plus two function keys for quick temperature setting and menu navigation
- User calibration facility for optimum accuracy at the required operating temperature
- Powerful integral pump – allows temperature-controlled fluid to be circulated to external equipment (16L/min, 210mbar)
- Dual-position bridge plate – ensures visibility/accessibility of the thermostat whilst optimising bench space
- Raised feet – for carrying / repositioning and retort stand access



- Liquid level protection and adjustable over temperature cut-out
- Clear 4 digit display – easy to read from a distance for instant reassurance
- Operating setpoint plus 3 adjustable temperature presets for convenience
- Robust construction, corrosion resistant materials, stainless steel tank – durable in demanding environments
- Excellent temperature stability and uniformity ensured by stirred circulation in the bath
- Drain tap allows easy emptying
- Choice of 120 V and 230 V models
- Optional insulated gabled, removable hinged lid designed to improve energy efficiency and prevent evaporation



\* see summary table on pp. 1.6–1.7 for accessories and for other models utilising the TC120 thermostat  
 \*\* operation below ambient temperature requires accessory cooling

- Applications:**
- Clinical, Microbiology and Pathology labs - media tempering, thawing & incubating samples
  - University research - temperature control of spectrophotometers & refractometers and jacketed vessels
  - Industrial labs - temperature probe calibration, water analysis, QC testing product, petrochemical testing, material testing, milk sample testing

## Heated circulating baths » TXF200-ST26 high specification showcase

## showcase 2 – high specification example

Model TXF200-ST26\* range -15°C to 200°C\*\*, stability  $\pm 0.01^\circ\text{C}$

High specification model with high performance digital thermostat and stainless steel tank for sophisticated applications requiring complex programming and/or ultra precise temperature control

- Optima™ high performance digital thermostat (TXF200) for ultra precise temperature control
- Stability  $\pm 0.01^\circ\text{C}$
- Cooling/heating range -15°C to 200°C\*\*
- Full colour screen
- Easy to program via interface or remotely via PC / Laptop using Labwise™ software
- Key functions easily accessed via home screen icons

TXF200-ST26 model shown

**Full colour screen** – clearly displaying actual and set temperatures, pump speed and clear status icons

**Intuitive screen icons and menus** – allow fast and accurate setup

**Socket for optional external probe** – allows remote temperature control

**Five-point user calibration facility** for optimum accuracy

**Countdown timer with audible alarm** alerts when your attention is required

**Drain tap** allows easy emptying

**Raised feet** – for carrying / repositioning and retort stand access.

**Memory capacity for 10 programs** containing 100 segments

**Program via intuitive user interface** or connect to PC/laptop to program via Labwise™ software

**The programming interface includes set target temperature** - a choice of time to target temperature or temperature ramp speed. An additional programmable relay for on/off control of ancillary equipment

**High and low temperature alarm settings** – visual, audible and programmable

**Powerful integral pump for external fluid circulation** – variable speed, 22L/min, 530mbar

**Optional insulated gabled and removable hinged lid** designed to improve energy efficiency and prevent evaporation



**Accessory cooling systems** allow operation at or below ambient temperature. See page 1.7 for details



\* see summary table on p. 1.6–1.7 for accessories and other models utilising the Grant high performance digital control units  
 \*\* operation below ambient temperature requires accessory cooling

#### Applications:

- Industrial labs - thermostat calibration, haze analysis (brewing), temperature probe calibration and material testing
- University research - temperature control of external equipment such as spectrophotometers and refractometers. Circulation of temperature control fluid to jacketed vessels

Heated circulating baths » T100-P12 budget showcase

## showcase 3 – budget example

Model T100-P12\* range ambient +5°C to 99°C, stability ±0.05°C

Economy model with digital thermostatic control unit and plastic tank for straightforward applications requiring accurate temperature control.

- **Optima™ digital thermostat (T100) for accurate temperature control**
- **Cooling/heating range ambient +5°C to 99°C**
- **Stability ±0.05°C**
- **3 programmable temperature presets**
- **Low liquid protection and fixed over temperature cut-out**



T100-P12 model shown

**Visual alarm** – alerts you when your attention is required

**Simple-to-use rotary dial and two function keys** for quick temperature setting and menu navigation

**Optional removable flat lid** to minimise evaporation of fluid and avoid contamination of samples

Choice of **120 V and 230 V models**

**User calibration facility** for optimum accuracy at the required operating temperature

**Low liquid protection and fixed over temperature cut-out**

Operating setpoint plus **3 adjustable preset temperatures** for convenience

**Wide range of optional accessories** for different applications, including test tube racks and raised shelves

**Robust plastic construction**, double-walled for rigidity, easy to clean

\* see summary table on p. 1.6 for accessories and for other models utilising T100 control units and/or plastic tanks

- Applications:**
- Clinical, Microbiology and Pathology labs - media tempering, thawing & incubating samples
  - Teaching labs, higher education/universities - practical demonstration/experimentation, sample preparation

## Heated circulating baths » Models, options and accessories

### Heated circulating baths – models, options and accessories

Any of the four Grant Optima™ digital thermostats can be combined with any of the Grant stainless steel and plastic tanks. The colour-coded summary table shows you the temperature range of each combination. For more details of Grant Optima™ thermostats see, p 1.8



Key to symbols	Heating circulators			
	General purpose digital		Advanced digital	
	T100	TC120	TX150	TXF200
<ul style="list-style-type: none"> <li> fixed over temperature cutout</li> <li> adjustable over temperature cutout</li> <li> display</li> <li> audible alarm</li> <li> timer</li> <li> pump</li> <li> external probe</li> <li> USB + RS232</li> <li> 2 2 point recalibration</li> <li> relay</li> <li> visual alarm</li> <li> 5 5 point recalibration</li> <li> menu system</li> <li> program storage</li> <li> programmable</li> </ul>	 <p>h: 335mm d: 172 mm w: 120 mm 2.5 kg</p>	 <p>h: 335 mm d: 172 mm w: 120 mm 2.5 kg</p>	 <p>h: 345 mm d: 172 mm w: 120 mm 3 kg</p>	 <p>h: 345 mm d: 172 mm w: 120 mm 3 kg</p>

### Tanks

Capacity (L) Outer tank dimensions	Working area (l x w) Min/max liquid depths Inner tank dimensions (l x w x h) Overall dimensions incl. controller (l x w x h)	2	2	5	
<b>ST5 – 5 L</b> stainless steel 3 kg h: 200 mm l: 330 mm w: 180 mm	<ul style="list-style-type: none"> <li>150 x 150 mm</li> <li>85/140 mm</li> <li>300 x 150 x 150 mm</li> <li>330 x 180 x 395 mm</li> </ul>	T100-ST5 amb.+15 to 100°C	TC120-ST5 0 to 120°C	TX150-ST5 0 to 150°C	TXF200-ST5 0 to 200°C
<b>ST12 – 12 L</b> stainless steel 4.5 kg h: 200 mm l: 360 mm w: 330 mm	<ul style="list-style-type: none"> <li>205 x 300 mm</li> <li>85/140 mm</li> <li>325 x 300 x 150 mm</li> <li>360 x 330 x 395 mm</li> </ul>	T100-ST12 0 to 100°C	TC120-ST12 0 to 120°C	TX150-ST12 0 to 150°C	TXF200-ST12 0 to 200°C
<b>ST18 – 18 L</b> stainless steel 7 kg h: 200 mm l: 540 mm w: 330 mm	<ul style="list-style-type: none"> <li>385 x 300 mm</li> <li>75/130** mm</li> <li>505 x 300 x 150 mm</li> <li>540 x 330 x 395 mm</li> </ul>	T100-ST18 0 to 100°C	TC120-ST18 0 to 120°C	TX150-ST18 0 to 150°C	TXF200-ST18 0 to 200°C
<b>ST26 – 26 L</b> stainless steel 7.5 kg h: 255 mm l: 540 mm w: 330 mm	<ul style="list-style-type: none"> <li>385 x 300 mm</li> <li>125/180** mm</li> <li>505 x 300 x 200 mm</li> <li>540 x 330 x 405 mm</li> </ul>	T100-ST26 0 to 100°C	TC120-ST26 -15 to 120°C	TX150-ST26 -15 to 150°C	TXF200-ST26 -15 to 200°C
<b>ST38 – 38 L</b> stainless steel 11 kg h: 255 mm l: 730 mm w: 330 mm	<ul style="list-style-type: none"> <li>575 x 300 mm</li> <li>125/180** mm</li> <li>690 x 300 x 200 mm</li> <li>730 x 333 x 405 mm</li> </ul>	T100-ST38 0 to 100°C	TC120-ST38 -15 to 120°C	TX150-ST38 -15 to 150°C	TXF200-ST38 -15 to 200°C
<b>P5 – 5 L</b> plastic 2.5 kg h: 180 mm l: 240 mm w: 330 mm	<ul style="list-style-type: none"> <li>120 x 150 mm</li> <li>85/140 mm</li> <li>240 x 160 x 150 mm</li> <li>390 x 200 x 380 mm</li> </ul>	T100-P5 amb.+15 to 99°C	TC120-P5 amb.+15 to 99°C	TX150-P5 amb.+15 to 99°C	TXF200-P5 amb.+15 to 99°C
<b>P12 – 12 L</b> plastic 3.5 kg h: 180 mm l: 415 mm w: 350 mm	<ul style="list-style-type: none"> <li>210 x 280 mm</li> <li>85/140 mm</li> <li>325 x 280 x 150 mm</li> <li>415 x 350 x 380 mm</li> </ul>	T100-P12 amb.+5 to 99°C	TC120-P12 amb.+5 to 99°C	TX150-P12 amb.+5 to 99°C	TXF200-P12 amb.+5 to 99°C
<b>P18 – 18 L</b> plastic 5 kg h: 180 mm l: 600 mm w: 365 mm	<ul style="list-style-type: none"> <li>280 x 325 mm</li> <li>85/140 mm</li> <li>510 x 290 x 150 mm</li> <li>600 x 350 x 380 mm</li> </ul>	T100-P18 amb.+5 to 99°C	TC120-P18 amb.+5 to 99°C	TX150-P18 amb.+5 to 99°C	TXF200-P18 amb.+5 to 99°C

Note: operation at or below ambient temperatures requires accessory cooling or a refrigeration unit on page 2.6




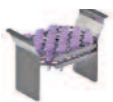

















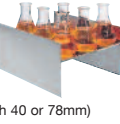


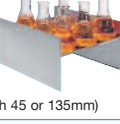













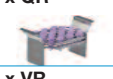

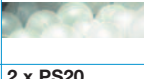






### Options and accessories

Labwise™ PC software (optional)				
Allows two-way communication for status display, programming and data capture (see p. 3.1 for more information) USB/RS232 cables provided	-	-		
<b>External probes</b> (optional) for monitoring and controlling temperature of remote loads				
<b>TXPEP</b> flexible plastic probe, 3m cable	-	-		
<b>TXSEP</b> stainless steel probe, 3m cable	-	-		
<b>Remote switching device</b> (optional)				
For switching appliances on and off (up to max. 8 Amps)	-	-	1	1
<b>Vertical turbine pumps</b> (optional)*				
Low noise, compact design. Supplied with pipe connections and special lid for fitting to tank, pipe bore 12.7 mm				
<b>VTP 1</b> max. pressure 1000 mbar max. flow 9 L/min  <b>VTP 2</b> max. pressure 1650 mbar max. flow 12 L/min	  	Required only where application demands a higher pressure than that delivered by the internal pump to maintain flow		

\* when pump is fitted, available working area is reduced \*\* maximum depth can be increased by 10 mm, by removing the circulation tray in 18, 26, 38 litre baths, with slight loss of performance

## Heated circulating baths » Options and accessories
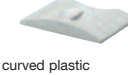
### Accessories

Lids*	Lids	Polypropylene spheres*	Rack systems†	Raised shelves	Accessory cooling systems**		
to help reduce evaporation/ heat loss and avoid sample contamination	For continuous use with water above 90°C. Stainless steel.	300 spheres in one pack (no. of packs required)	to optimise use of available bath capacity (no. of racks accommodated)	to allow shallow vessels to be accommodated	to allow systems to operate at or below ambient temperature by means of a cooling coil dipped into the bath; designed for minimal impact on working area		
					<b>Refrigerated immersion coolers</b> Consist of a cooling coil connected to a refrigeration unit by a flexible pipe. Extract heat continuously, with the bath control unit controlling temperature 	<b>Heat exchange coil</b> Designed to be attached to a supply of cooling tap water or a refrigerated circulator	
					<b>C1G</b> (0 to 40°C***)	<b>C2G</b> (- 15 to 40°C***)	<b>CW5</b> (2°C above coolant temperature)
<b>ST5</b>  flat stainless steel	-	1 x PS20 	1 x QR 	-		-	
<b>ST12</b>  gabled, hinged (removable) stainless steel	LST12 	1 x PS20 	2 x VR 	RS14  (h 40 or 78mm)		-	
<b>ST18</b>  gabled, hinged (removable) stainless steel	LST26 	2 x PS20 	4 x VR 	RS22  (h 40 or 78mm)		-	
<b>ST26</b>  gabled, hinged (removable) stainless steel	LST26 	2 x PS20 	4 x VR 	RS28  (h 45 or 135mm)			
<b>ST38</b>  gabled, hinged (removable) stainless steel	LST38 	3 x PS20 	6 x VR 	RS28 or RS38  (h 45 or 135mm)			
<b>P5</b>  flat, stainless steel	-	1 x PS20 	1 x QR 	-	-	-	-
<b>P12</b>  curved plastic	-	1 x PS20 	2 x VR 	RS14  (h 40 or 78mm)	-	-	-
<b>P18</b>  curved plastic	-	2 x PS20 	4 x VR 	RS22  (h 40 or 78mm)	-	-	-

\* Between operating temperatures 60°C and 100°C and below room temperature a lid or layers of polypropylene spheres should be used.

\*\* The cooling coil can be continuously immersed in liquids up to 100°C with the cooler switched off, and may be used to cool liquid down from 100°C, but it is not designed for continuous operation above 40°C.

\*\*\* Minimum operating temperature without accessory cooling is ambient + 5°C (amb.+ 15°C for P5 and ST5 tanks).

† Rack capacity (no. of test tubes per rack)

VR racks	Tube size	Capacity	QR racks	Tube size	Capacity
VR-13	ø 10-13 mm	65	QR-13	ø 10-13 mm	30
VR-19	ø 16-19 mm	36	QR-19	ø 16-19 mm	16
VR-24	ø 24 mm	23	QR-24	ø 24 mm	10
VR-30	ø 30 mm	14	QR-30	ø 30 mm	5
VR-SE	0.5 ml	102	QR-SE	0.5 ml	44
VR-LE	1.5 ml	75	QR-LE	1.5 ml	35









## Heated circulating baths » Technical specifications

## Heated circulating baths – technical specifications

## Grant Optima™ thermostats

● = standard

			General purpose digital		Advanced digital	
			T100	TC120	TX150	TXF200
						
Stability (DIN 12876)@70°C	°C		±0.05	±0.05	±0.01	±0.01
Uniformity (DIN 12876)@ 70°C	°C		±0.1	±0.1	±0.05	±0.05
Setting resolution	°C		0.1	0.1	0.1 (0.01 with Labwise)	
Display			4 digit LED		full colour QVGA TFT	
Timer function			–	1 to 6000 mins	1 min to 99 hrs 59 mins	
No. of temperature presets			3	3	3	3
Re-calibration points			2	2	5	5
Socket for external probe (TXPEP, TXSEP)			–	–	●	●
Communications interface			–	–	USB, RS232	USB, RS232
Programmable			–	–	remote via PC / laptop 1 program / 30 segments	direct via user interface or remote via PC / laptop 10 programs / 100 segments
Relays			–	–	1	1
Safety	over temperature		fixed		adjustable cut-out	
	fluid level – float		●	●	●	●
Language capability			–	–	EN, FR, DE, IT, SP	EN, FR, DE, IT, SP
Alarms (can be configured to switch a relay)			–	high (no relay)	high and low	high and low
Heater power	230 V	kW	1.3	1.3	1.9	1.9
	120 V	kW	1.4	1.4	1.4	1.4
Electrical power	230 V	kW	1.4 (50-60 Hz)	1.4 (50-60 Hz)	2.0 (50-60 Hz)	2.0 (50-60 Hz)
	120 V	kW	1.5 (50-60 Hz)	1.5 (50-60 Hz)	1.5 (50-60 Hz)	1.5 (50-60 Hz)
Height above tank rim		mm	200	200	200	200
Depth below tank rim		mm	135	135	145	145

## Grant Optima™ thermostat pumps (integral)

Maximum pressure	water	mbar	–	210	310	530
Maximum flow	water	L/min	–	16	18	22 (adjustable flow rate)
Pump connector	6 mm bore*		–		fits 9 mm inner diameter tubing	
Pump connector	11 mm bore*		–		fits 15 mm inner diameter tubing	







\* 6 and 11 mm bore pump connectors supplied as standard. For more options see page 1.9

Grant immersion thermostats are suitable for use with Grant stainless steel and plastic tanks. With the addition of a clamp (T-clamp) they can also be attached to virtually any vertical sided tank with a maximum wall thickness of 35 mm for rectangular tanks, 30mm for circular tanks (300 mm diameter), and a capacity of up to 50 litres. Minimum and maximum temperatures achievable are dependent upon the tank insulation and minimum operating temperature depends on the accessory cooling device.






## Heated circulating baths » Technical specifications

### High pressure pumps (optional)

			VTP pumps						
			VTP1-FLST5	VTP1-FLST12	VTP1-FLST26	VTP2-FLST5	VTP2-FLST12	VTP2-FLST26	
									
Fits ST tank			ST5	ST12	ST18/26	ST5	ST12	ST18/26	
Maximum pressure	water	mbar	1000			1650			
Maximum flow	water	L/min	9			12			
Pipe bore	inlet/outlet	mm	12.7			12.7			
Electrical connection	10 amp IEC			10 amp IEC					
Power consumption		W	30			40			
Power output to liquid @ 20°C		W	15*			22*			
Safety	thermal fuse			thermal fuse					

### Accessory cooling systems

			Immersion coolers		Heat exchange coil
			C1G	C2G	CW5
					
Cooling power	@ 20°C	W	350	400	-
	@ 0°C	W	110	320	-
	@ -10°C	W	-	170	-
Overall consumption		VA	300	500	-
Dimensions	d/w/h	mm	460/305/225		-
Weight		kg	17	21	0.1
Flexible pipe	l	mm	925	925	-
Coil	ø / l	mm	77/55	77/55	77/55
Pipe bore inlet/outlet		mm	-	-	7
Electrical supply	120 V (60 Hz) or 230 V (50Hz)				

\* The VTP optional pumps will transfer additional heat to the baths, so the minimum temperature achievable with or without accessory cooling will be increased. Note: when ordering a VTP pump, please specify which Grant tank it is to be used with.

### Pump connectors (optional)

	Part number
Replacement plastic pump inlet/outlet connector. Fits tubing 9mm inner dia. Temperature range -50 to 200°C	P-M6
Replacement plastic pump inlet/outlet connector. Fits tubing 15mm inner dia. Temp range -50 to 200°C	P-M11
Stainless steel pump inlet/outlet connector, M16 x 1 male. Fits M16 hose. Temp range -50 to 200°C	M-M16
Metal pump inlet/outlet connector, dual seal super rapid 4mm. Fits semi rigid tubing 4mm outer dia. Temp range -20 to 100°C	M-SR4
Metal pump inlet/outlet connector, dual seal super rapid 6mm. Fits semi rigid tubing 6mm outer dia. Temp range -20 to 100°C	M-SR6
Metal pump inlet/outlet connector, dual seal super rapid 8mm. Fits semi rigid tubing 8mm outer dia. Temp range -20 to 100°C	M-SR8
Metal pump inlet/outlet connector, hose barb 7mm. Fits flexible tubing 7mm inner dia. Temp range -40 to 120°C	M-HB7
Metal pump inlet/outlet connector, hose barb 9mm. Fits flexible tubing 9mm inner dia. Temp range -40 to 120°C	M-HB9
Metal pump inlet/outlet connector, hose barb 12mm. Fits flexible tubing 12mm inner dia. Temp range -40 to 120°C	M-HB12
Metal pump inlet/outlet plate, 1/4 " BSP/G1/4 female. Temp range -50 to 200°C	M-UC

# High temperature oil baths / circulators

## HED series


Series of robust baths/circulators for high temperature applications combining electronic control with Grant's proven circulation system to provide long-term stable, uniform temperature control. Designed for use with high temperature liquids.

- Temperature range 50°C to 260°C
- Stability  $\pm 0.25^\circ\text{C}$ , uniformity  $\pm 0.2^\circ\text{C}$
- Timed or continuous operation
- Choice of two tank sizes – 10 litres (HE10D) and 30 litres (HE30D)

**Insulated flat lid** to ensure user safety. Model HE30D has a 2-part lid to facilitate easy access

**Optional accessory pump HQ** available for external circulation (closed system only) – mounted on lid for convenient fitting and use

**Robust construction for durability and heat efficiency** – heavily insulated stainless steel tank in a stove enamelled outer case. Exposed or immersed parts are made from corrosion resistant materials

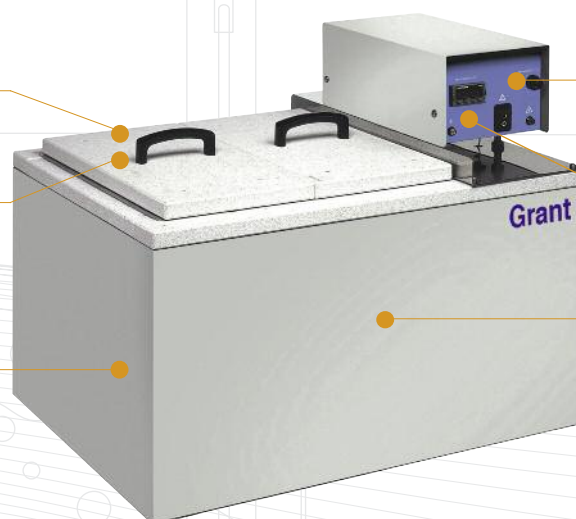

**Digital controller raised** to allow for cooling airflow above the tank

**Fitted with extra powerful stirrer and a heater** with a large surface area to reduce local overheating and deterioration of the oil

**Inbuilt safety features** ensure that the set temperature is not exceeded and the bath does not operate without a minimal level of liquid



### Accessory pump for external circulation (closed systems only)

- **HQ10** – for 10 litre bath HE10D
- **HQ30** – for 30 litre bath HE30D

- Applications:**
- Academia - crude oil extraction
  - Industry - oil and gas research

Heated circulating baths » HED series » Models and specifications

Products for special high temperature applications – models and specifications

			High temperature baths/circulators digital control	
			HE10D	HE30D
				
Temperature range	°C		50 to 260	
Stability (DIN 58966)	°C		±0.25 (at 150°C)*	
Uniformity	°C		±0.2 (at 150°C)*	
Display			LED	
Display resolution	°C		1	
Timer	mins		1 to 9999	
Alarms			high/low	
Heater power 230 V	kW		1.1	2.2
Overall consumption 230 V	kW		1.2	2.3
Liquid flow rate, maximum	L/min		6**	
Pump head pressure @ 0 L/min	m		1.6**	
Pipe bore	mm		7.5*	
Working area/tank opening	mm		190 x 210	350 x 305
Tank dimensions l x w x h	mm		305 x 210 x 185	455 x 305 x 240
Liquid depth min/max	mm		100/140	145/190
Safety	over temperature protection		adjustable cut-out	
	liquid level		float switch	
Electrical power	230 V 50/60 Hz	kW	1.2	1.3
	120 V 50/60 Hz	kW	1.2	1.6
EMC emissions			Class A	
Weight	kg		19	30

\* using silicon oil # using accessory pump

## Heated circulating baths » Gerber bath

# Gerber Bath

Robust, high precision temperature-controlled stirred bath specifically designed for use in dairy laboratories, with sufficient depth to allow immersion of Gerber butyrometers. Also suitable for general laboratory applications where a deeper tank is required.

- Temperature range ambient +5°C to 100°C
- Stability  $\pm 0.05^\circ\text{C}$
- Precision digital thermostatic temperature control, with stirred circulation
- Capacity for up to 36 Gerber butyrometers
- Fixed over temperature cutout
- 3 programmable temperature presets

**Removable control unit** for easy cleaning and servicing. All immersed parts – the heater, float, stirrer and temperature sensor – are made of materials that will resist occasional contamination of the water by spilt Gerber acid

**Space for three purpose-designed racks** (H1-25M), each with the capacity to hold 12 butyrometers

**Robust stainless steel inner tank and outer case** with a hinged lid plus handle for convenient access

**User calibration facility** for optimum accuracy at the required operating temperature



## T100-MST20 Gerber bath

**The bright LED display** constantly displays actual temperature for at-a-glance reassurance

**Sophisticated digital control unit** provides adaptive intelligent PID temperature control

**Low liquid level protection and fixed over temperature protection**

**Simple-to-use rotary dial plus two function keys** for quick temperature setting and menu navigation

**Applications:** ■ Gerber testing to determine the fat content in milk and dairy products

Heated circulating baths » Gerber bath » Models and specifications

Gerber bath – models and specifications

Gerber bath

T100-MST20



h: 380 mm  
d: 600 mm  
w: 1040 mm

Temperature range	°C	ambient +5 to 100	
Stability	°C	±0.05	
Display		4-digit LED	
Display resolution	°C	0.1	
Heater power	230 V	kW	1.3
	120 V	kW	1.4
Working area/tank opening	mm	240 x 210	
Tank capacity (nominal)	L	20	
Liquid depth	min/max	mm	185/225
Electrical power	230 V 50/60 Hz	kW	1.4
	120 V 50/60 Hz	kW	1.5

