
Important Safety Information

- Read the full manual before use. Safety may be impaired if it is not followed.
- Only connect to the voltage and frequency shown on the serial number plate. An earth is required.
- Always ensure the bath is filled with a liquid rated for the working temperature. See manual for liquid selection.
- Ensure the TX150/TXF200 is positioned so that the plug is easily accessible for disconnection.
- Before moving or cleaning, allow the unit to cool down and disconnect from the power supply.
- Safety can only be assured if the TX150/TXF200 is securely installed on a suitable vessel. See manual for guidance.

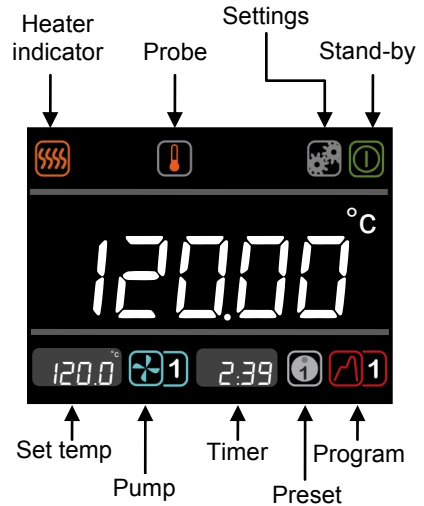Setting the temperature


Rotate the dial until the set temperature icon is highlighted.
 Press the **S** button to select. Rotate the dial to the desired temperature.
 Press **S** to set.

Setting the over-temp

Rotate dial to desired temp cut-out (above set point)

Using the external pump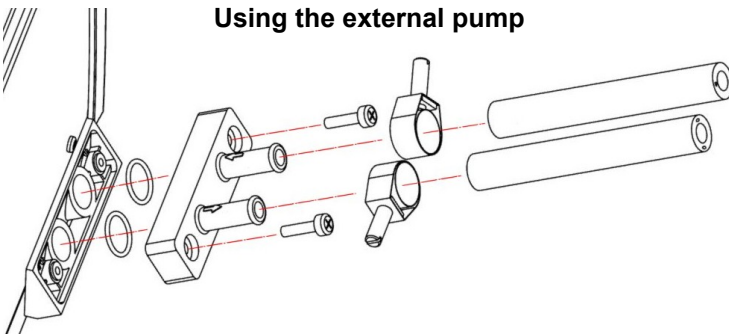

Ensure the TX150/TXF200 is switched off. Connect hoses securely to the pump outlet plate as shown left.

Always use pump connectors and hoses that are suitable for the operating temperature and liquid used.

Never disconnect any pipes or hoses while they contain very hot or very cold liquids or while the unit is pumping.

Safety cut out and alarm light

The TX150/TXF200 is fitted with a float switch which monitors the liquid level during operation. Should the liquid level drop below a certain level, the safety cut out will be activated, and alarm light will appear. Top up the bath with liquid until at least a minimum level to resume operation.

Cleaning

Clean the unit with a damp cloth, using only water. Do not use chemical cleaning agents.

Consult manual for full instructions.

Service & repair

For service or repair please contact our Service Department or your local distributor. All returned units must be accompanied by a Return Materials Authorisation (RMA) number, obtainable by contacting the Grant service department.

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